SONICA SOUNDS BETTER.

NEW DESIGN
NEW TECHNOLOGY
NEW LIFE.

SOLTEC website

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SOUNDS BETTER.

ITS PRODUCTS

SOLTEC designs and produces ultrasonic cleaners and detergents for professional use in all fields in which high cleaning standards must be achieved within the shortest possible timeframe such as in the chemical, electronic, pharmaceutical, dental, medical, military and aerospace sectors. Using ultrasonic energy and a process known as cavitation, ultrasonic cleaning is the most effective way of deep cleaning parts that are difficult to reach as well as complex shaped items (mechanical parts, gears, moulds, precision instruments).

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SOUNDS BETTER.

Ultrasonic cleaners

At present **SOLTEC** produces over 20 models with more than 50 different versions.

Each component was designed and produced for achieving optimal results in the field of precision cleaning. Used in many sectors, both at an industrial level and in individual laboratories, **SONICA** ultrasonic cleaners are specifically made to order and can be designed in a standard version or special version that complies with the specific requirements and needs of individual customers.

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1. Piezoelectric Transducers.
2. Sweep System oscillator module.
3. Microprocessor circuit base for degassing, heating and ultrasonics (EP Version).
4. Stainless steel nipple TIG welded providing a watertight seal.
5. Stainless steel with special surface treatment.
6. Waterproof membrane keyboard with led display (EP Version).
7. Drain Cock 1/2".
8. Sturdy handles.

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The new **SONICA S3** range of ultrasonic cleaners is characterized by its newly-designed control panels. The innovative layout has been designed to be user friendly.

All **SONICA S3** cleaners with a capacity of 3 litres or more have been equipped with state-of-the-art high-power piezoelectric transducers and with ultrasonic generators using the innovative SweepSystem oscillation technology, providing for perfect cavitation of detergents.

All new **SONICA S3** cleaners come with a 24 month warranty.

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On the left: SONICA 3300EP during ultrasonic cleaning.
Below: SONICA 4300 flush-mountable version.

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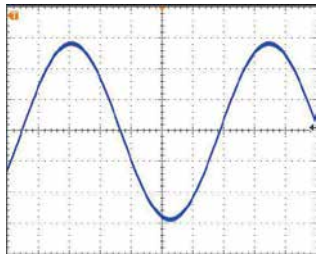
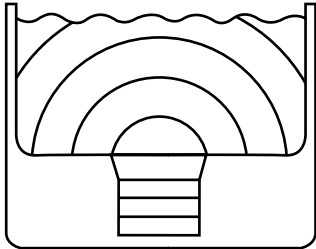
Comparison of acoustic field distribution inside the liquid.

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Comparison of the generators' electric output signals.

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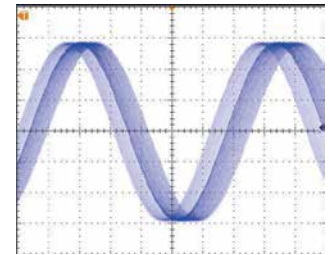
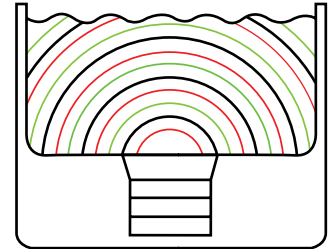
Comparison of irradiation uniformity using the two methods on a sample aluminium sheet.

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The new **SONICA S3** series with Sweep System technology offers the best performance available from ultrasonic cleaning equipment.

The state-of-the art generator incorporates a special ultrasonic oscillator using Sweep System Technology.

The frequency from the ultrasonic generator operates around a central frequency of 39 Khz and the transducers shift between 38 and 40 Khz. Sweep System Technology gives:

- Faster cleaning.
- Damage free cleaning.
- Increased distribution of ultrasound energy.
- Better cleaning.
- Improved cavitation.

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Ultrasonic precision cleaning is the fastest and most economical way of effectively removing oil, grease, filings, small chips, surface dust, biofilm, organic matter, etc. In the following fields of applications:

1. Mechanical Industry & Degreasing.
2. Electronic Industry.
3. Chemical Labs.
4. Cleaning of Metallographic, Ceramic, and Mineralogical Specimens.
5. Firearms Maintenance.
6. Jewels and Clock Industry.
7. Winch Maintenance.
8. SCUBA & SCBA cleaning.
9. Dental/Medical and Infection Control.

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Mod. M

The simplest version in the SONICA range is equipped with a timer which can be set for up to 15 minutes. It can also be set to manual operation. **Sweep System** generator included.

Mod. MH

This is equipped with a timer which can be set for up to 15 minutes and can also be set to manual operation. The thermostat for the heater is set at 60°C. **Sweep System** generator included.

Mod. ETH

Cleaning times and temperatures controlled electronically by a microprocessor. Waterproof control panel. **Sweep System** generator included.

Mod. EP

The most technologically advanced of the SONICA range: modern digital display control panel; degassing time setting; cleaning times from 0 to 99 minutes; tank liquid temperature setting up to 70°C; storage capacity for up to 10 cleaning programmes; **Jet program** function allows for quick and easy cleaning cycles. **Sweep System** generator included.

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 The **SONICA S3** industrial series includes ultrasonic cleaners with a tank capacity of 45 to 130 litres. Besides offering its standard models, SOLTEC can also design "Custom" ultrasonic cleaners, based on its customers' specific needs and requirements.

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* Available also in "i" integrated version.

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Technical specifications subject to change.

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Accessoires

1. Stainless steel holder basket.
2. Special stainless steel holder for volumetric flask.
3. Auxiliary stainless steel tank for indirect cleaning with aggressive detergent. Can also be used as a rinse or holding tank after the ultrasonic cleaning.
4. Small stainless steel immersion basket for cleaning small objects.
5. Glass beaker, complete with retaining ring.
6. Stainless steel lid.
7. Stainless steel cassettes carrier for dental instruments. Holds up 5 instruments.
8. Glass beaker positioner made in electro finished stainless steel.
9. SONICA decontamination tank kit: 1 stainless steel tank autoclavable; 1 stainless steel autoclavable lid; 1 metal positioning case for stainless steel tank; 1 SONICA CL 4% concentrated disinfecting detergent.

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The only automatic system operated by ultrasonics
and a Thermal Disinfection cycle at 93°C for the treatment of surgical instruments

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THE REAL INNOVATIVE SAFETY SYSTEM FOR PATIENTS AND OPERATORS ALL IN ONE APPLIANCE.

Ultrasonic cleaning with automatic detergent / disinfectant liquid dosage, multiple rinsing cycles with water supply, Thermodry treatment at 85°C, special Thermal Disinfection cycle at 93°C for 3 minutes, drying cycle and USB recording of the relative data of the surgical instruments treatment cycles **in an only appliance.**

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Ventilation system included in the cover

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Detergent/disinfectant liquid loading

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Liquid drain tube connected to an internal drain pump

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Spray nozzles for multiple rinsing cycles

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Innovative Thermodry system at 85°C
Thermal Disinfection cycle at 93°C for 3 minutes
Automatic multiple rinsing cycles (patents) Nr. CO2011A000008
Nr. CO2011A000009

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SONICA® TRENDMATIC innovative system is controlled by a 32 bit microprocessor and each phase of the instruments treatment can be set up by the final user through a wide LCD graphic touch-screen display.

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All rights reserved SOLTEC Srl
Variations may be made without notice.

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* Special detergent additive for Winch Cleaning 2 product - add 1% on Winch Cleaning 2 total amount - If it is used only as simple detergent dilute it with water from 5% to 10% depending on the extent of the deposits.

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SOUNDS BETTER.

 SONICA® Detergents

SONICA detergents have been specially designed to allow ultrasound waves to be propagated and ensure thorough and accurate cleaning and disinfecting. The range includes detergents and disinfectants that are able to remove all kinds of deposits from any type of surface. Directly formulated in the **SOLTEC** laboratories, detergents in the **SONICA** range can also be created using customized formulas that remove specific types of contaminants. All **SONICA** detergents are for strictly professional use and are subjected to laboratory tests. Guaranteed safe to use on treated surfaces and materials.

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New **Desafío Español 2007** team's official supplier at **32nd America's Cup Challenger**

SOLTEC, leading manufacturer of ultrasonic cleaning equipment **SONICA**, has become the Official Supplier of the America's Cup team Desafío Español 2007.

The Managing Director of the Desafío Español 2007 team, Mr. Agustín Zulueta and the Managing Director of **SOLTEC** Srl, Mr. Stefano Guidi, on 5th June 2006, signed an agreement in which **SOLTEC** Srl was appointed as 'Official Supplier' of the Desafío Español 2007 team.

Not only did **SOLTEC** provide the hardware for the process of winch cleaning, but **SOLTEC** also offered its technical background and expertise in ultrasonics with an 'on-site' technical team at the Desafío Español base. This was **SOLTEC**'s second experience at the America's Cup, during the 31st edition **SOLTEC** was OneWorld Challenge Team's Official Supplier.

SOLTEC supplied an 18lts capacity **SONICA** 4300EP featuring the most advanced technology available in ultrasonic tanks - 'Sweep System Generator', the winch team was able to clean components to a vastly improved level.

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MISSION. Its mission is to design and produce products and services in the field of precision cleaning based on its end customers' specific requirements and guarantee excellent quality/price ratio.

VISION. Its vision is to continue to improve its offer by achieving a high level of quality and low costs and low environmental impact that meet the demands and expectation of its customers and anyone else involved.

STRATEGY. To achieve these key objectives, **SOLTEC** focuses on drawing from the full potential of the managerial and technical expertise of its human resources and on creating a professional line of products using state-of-the-art technology.

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**Copyright © 1999 - 2017 Soltec S.r.l. tutti i diritti riservati.
SOLTEC S.r.l.